


Discussion


Identification and elimination of interface contaminants is critical for successful production and the ultimate performance of semiconductor devices. Interface contaminants include, but are not limited to, organic and inorganic cleaning residues, residual etch stops, and metals, and deposition chamber contaminants. In order to detect interface contamination on patterned devices an analytical technique must have good spatial resolution and good sensitivity for a wide range of elements. Numerous techniques are suitable for devices with features greater than $\sim 50\mu\text{m} \times 50\mu\text{m}$. For features smaller than $1\mu\text{m}$, FE-Auger is the technique of choice for analyzing interfacial contaminants.

The first example is a laser device that did not meet performance specifications. The facet coating was Au/Pt/Ti on GaAs. The region of interest is approximately 50 microns wide and several hundred microns long. An Auger depth profile was acquired from the center of the device as shown in the accompanying secondary electron image. Silicon was detected between the titanium layer and the GaAs substrate. The source of the silicon was identified as remnant of silicon nitride that was not completely removed prior to the deposition of the titanium. The profile also shows some oxygen present in the titanium layer near the GaAs interface.

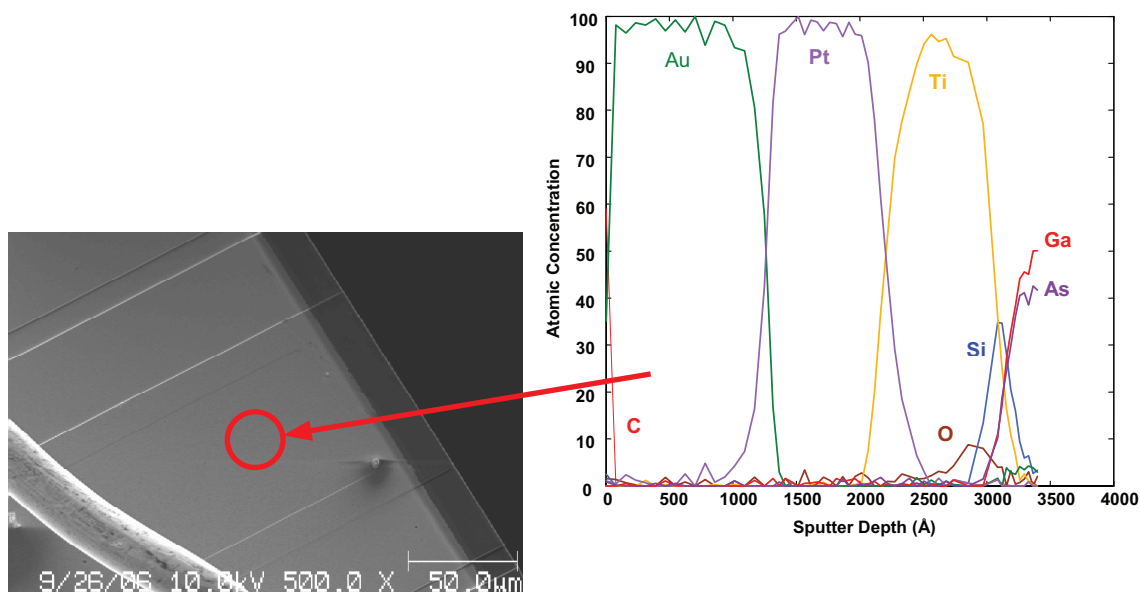


Example #1 Auger sputter depth profile acquired from a laser semiconductor device shows the presence of silicon at the Ti:GaAs interface. The source of the silicon was residual Si_3N_4 .

In the second example, the objective of the analysis was to identify contaminants at a silicon nitride-gallium arsenide interface within a $1\mu\text{m}^2$ area. The depth profile was stopped at the silicon nitride-gallium arsenide interface so that a survey spectrum could be acquired. The profile and interface survey spectrum revealed the presence of indium and phosphorous at the interface. The contamination was identified as residual InGaP which used as an etch stop layer that was not completely removed during wafer processing.

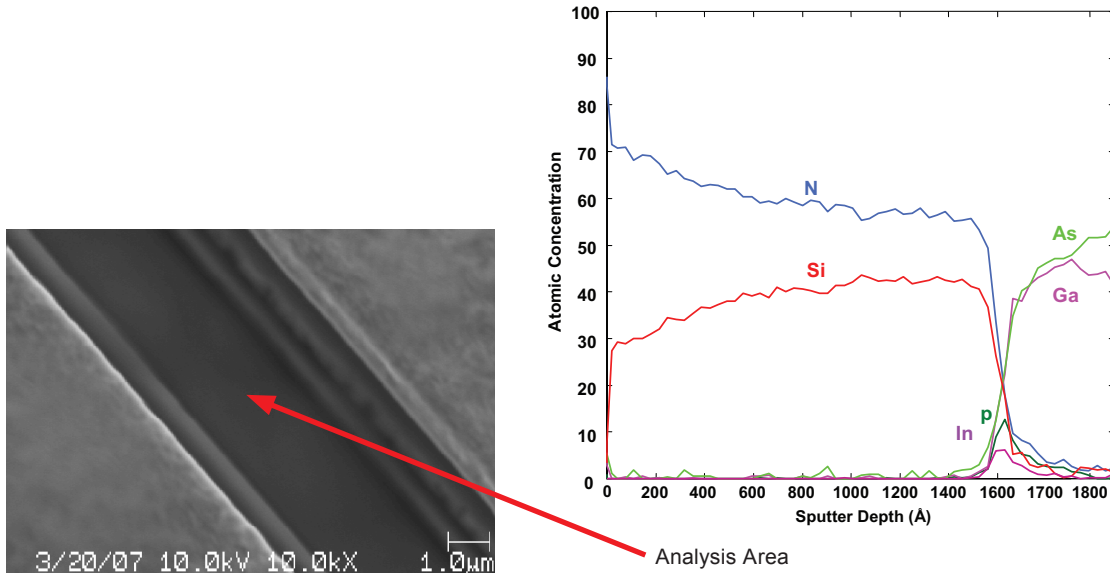

Example #2 Auger sputter depth profile acquired from a $1\mu\text{m}^2$ area shows the presence of indium and phosphorous at the $\text{Si}_3\text{N}_4:\text{GaAs}$ interface. The contamination was identified as residual InGaP etch stop.

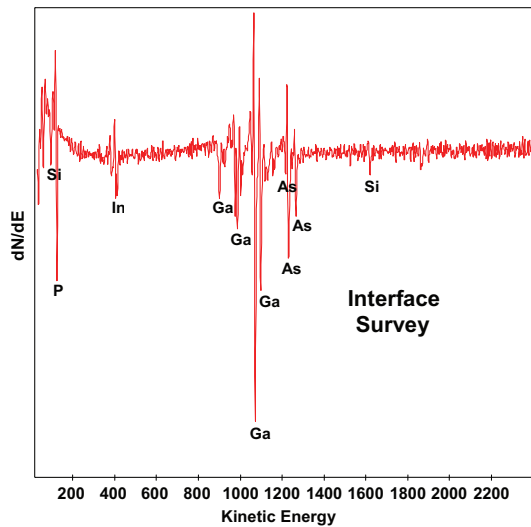

Visit www.eaglabs.com for more information about all of EAG's services and solutions.

EAG Limited Administrative Offices, 810 Kifer Road, Sunnyvale, CA 94086 phone: 408 530 3500

Copyright © 2008 EAG Limited. All rights reserved. EAG, the EAG logo, are registered trademarks of EAG Limited.

Evans Analytical Group, Charles Evans & Associates, Thin Film analysis, Inc., Applied Microanalysis Labs, Inc., AMIA Labs, Advanced Materials Engineering Research, Cascade Scientific Ltd., Cascade Scientific GmbH, Nano Science Corporation, Shiva Technologies, Inc., Shiva Technologies Europe SARL, Accurel Systems International Corporation, Micro Electronic Failure Analysis Services, Inc., DSL Labs Inc., White Mountain Labs LLC, are service marks of EAG Limited. All other company, product and service names may be trademarks of their respective companies. While every effort is made to ensure the information given is accurate, EAG Limited does not accept liability for any errors or mistakes which may arise. All information is subject to change without notice.